

Keys to AAV Manufacturing Success: Fast, Robust and High-yield Process Development with DoE, Automated Bioreactors and Selected Critical Raw Materials

Building a scalable, data-driven upstream platform for industrial AAV manufacturing

Hakima FLICI Scientific Support Specialist


Agenda

1. rAAV gene therapy: background and current challenges
2. Sartorius solutions for AAV process development
3. Case Study: Development of a robust rAAV upstream platform

Optimization of Bioreactor and transfection parameters (small scale)

Process scale-up


4. Key results and takeaways

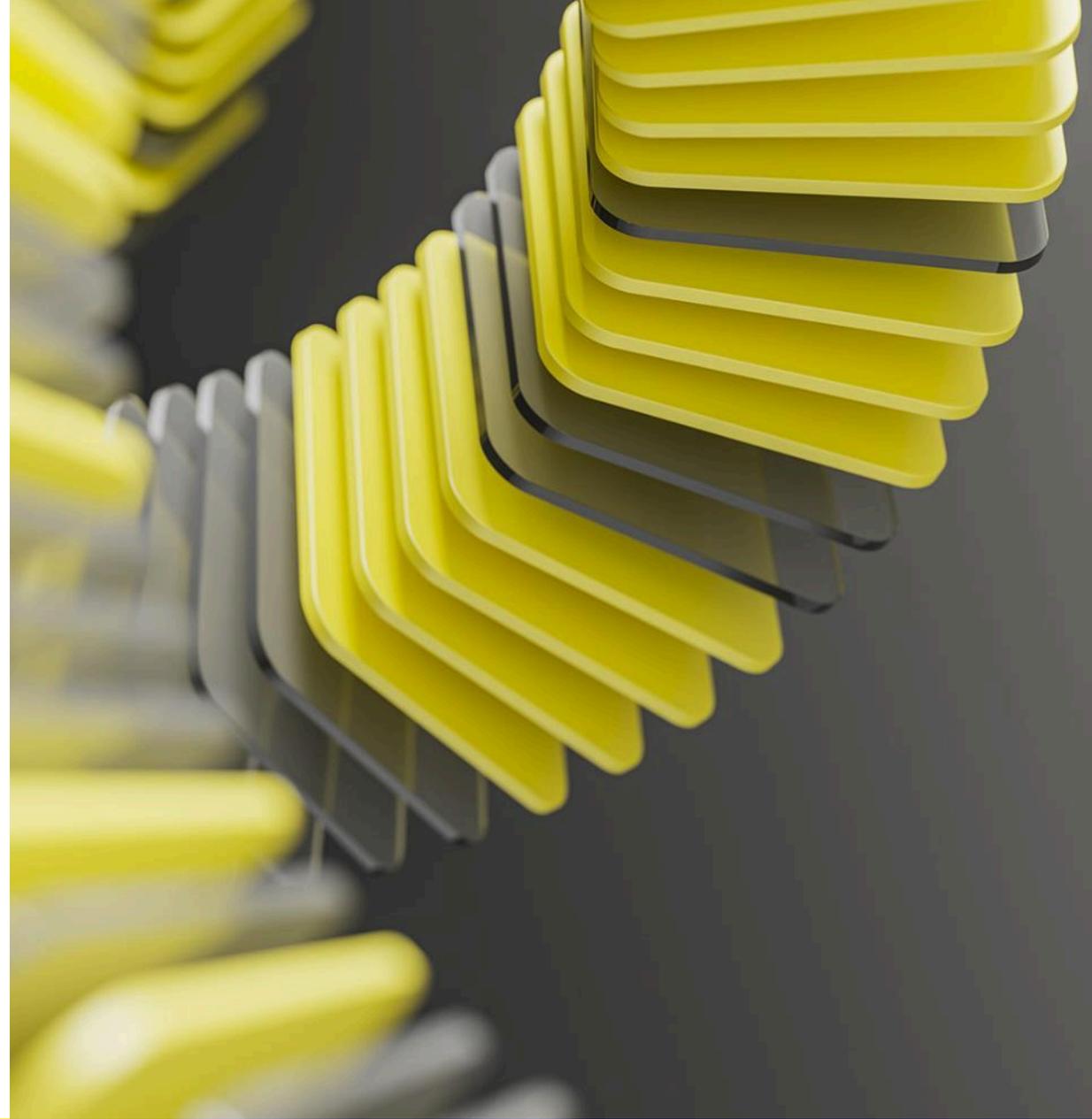


rAAV gene therapy: background and current challenges

Expanding Demand Highlights Urgency in AAV Gene Therapy Advancement


Unique AAV-based Drugs by Highest Development Stage

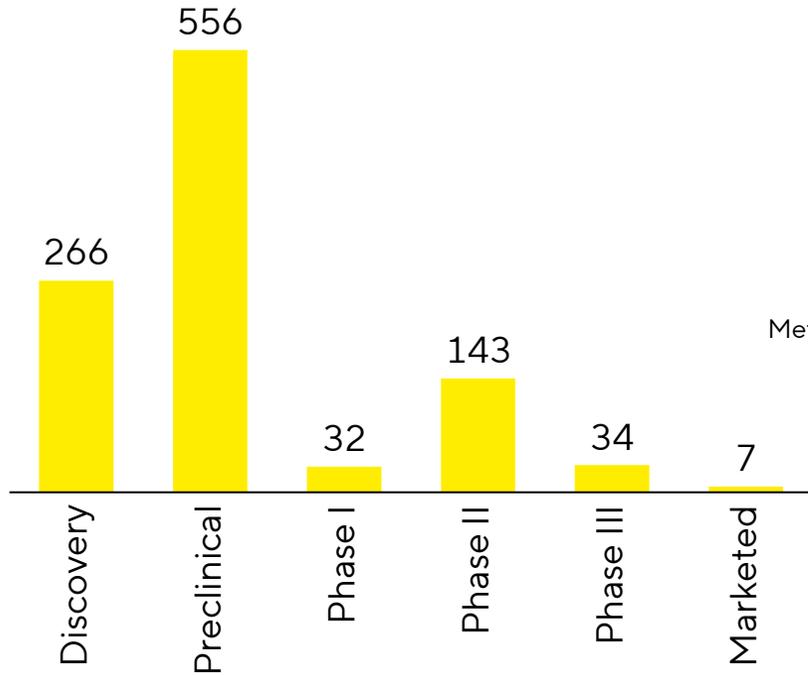


Unique AAV-based Drugs by Therapy Area

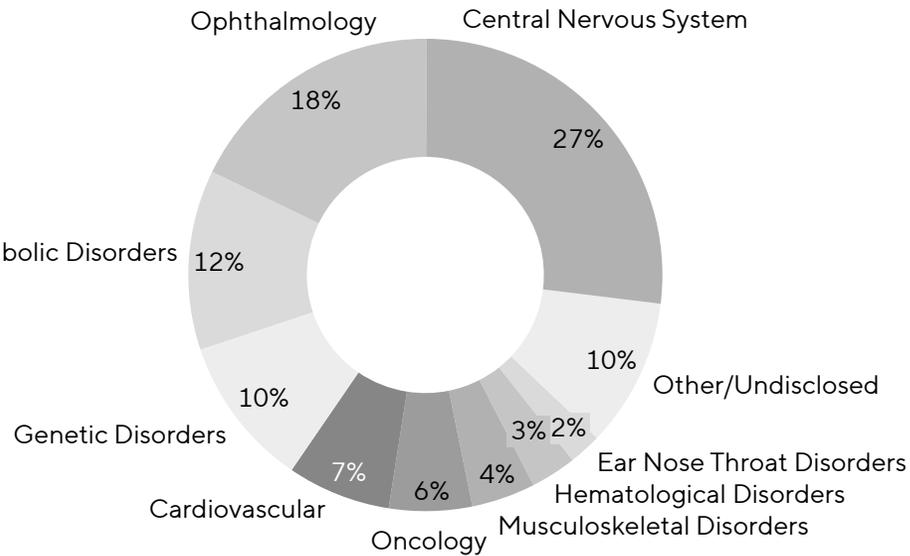

Proportion of different serotypes used

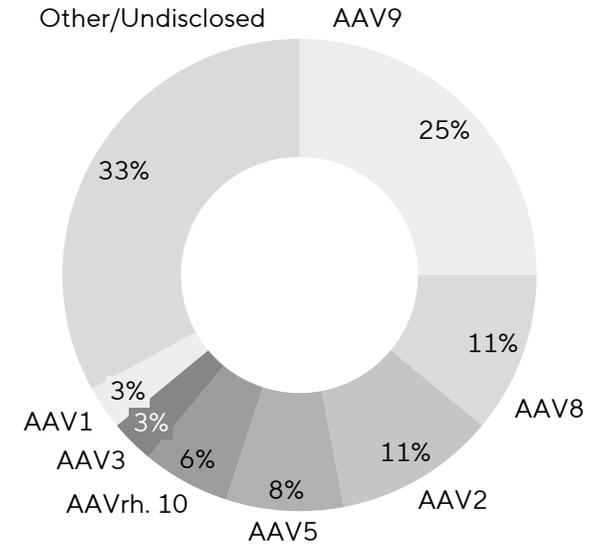

Overview of AAV Production Using Plasmid Transfection

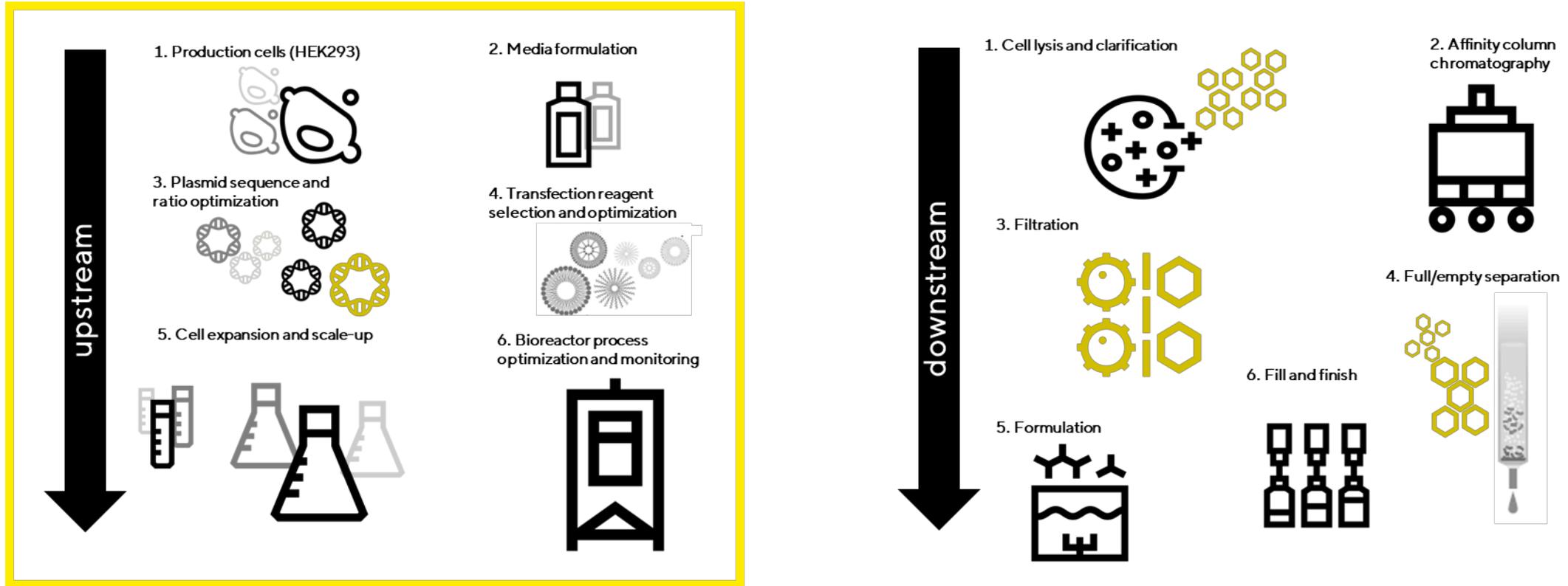


Adapted from *Cell & Gene Therapy Insights* 2024; 10(6), 821-840 DOI: 10.18609/CGTI.2024.095

Sartorius solutions for AAV process development

Advanced Therapies 2026 | Do not replicate

Comprehensive solutions for CGT development & manufacturing

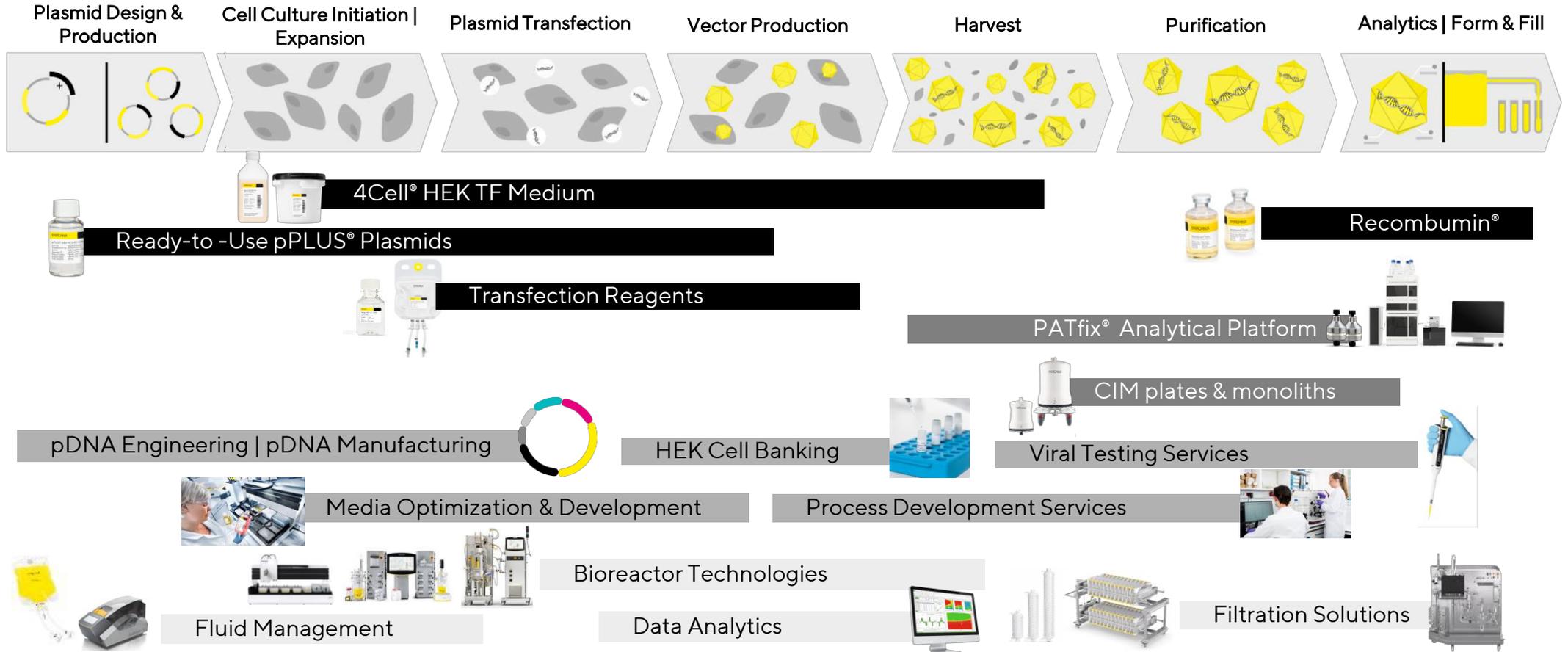

Case Study: Development of a robust rAAV upstream platform


Building a Robust AAV Upstream Platform

Objectives

Increase yield & Develop a Robust AAV Production Platform by Leveraging Comprehensive Design-of-Experiment (DoE) & High-Throughput Automated Bioreactors Powered by Next Gen Raw Materials




Raw Materials Used in This Study

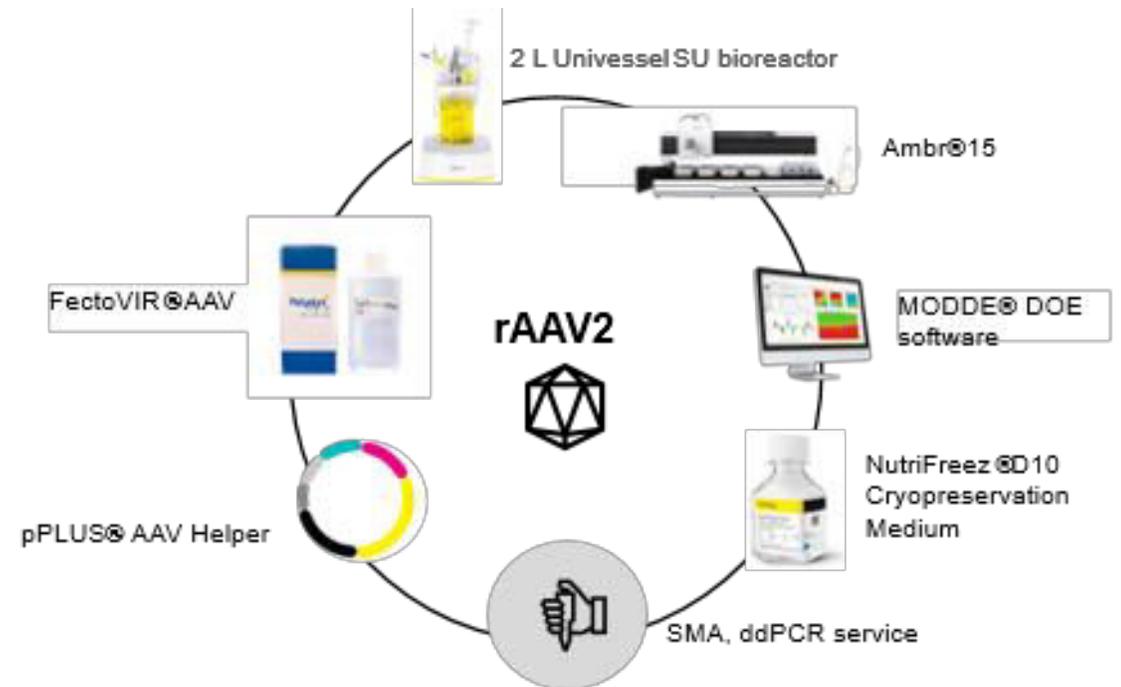

AAV Production, Where We Started

rAAV2-GFP : Productivity of $1.1\text{E}+10$ vg/mL (Bulk harvest) with the standard culture medium in shake flask.

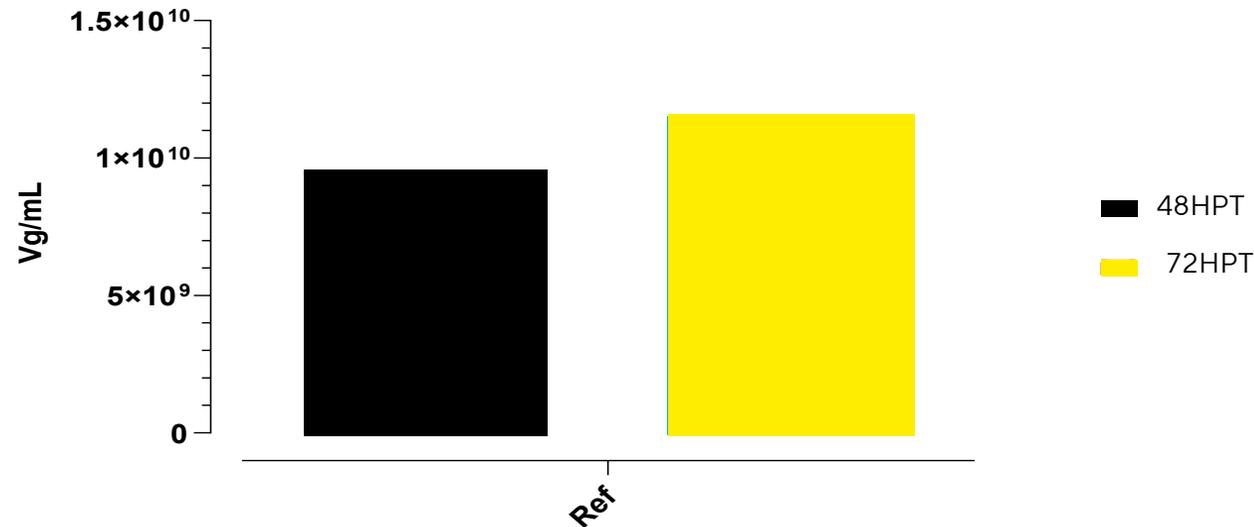

Significant productivity gains are needed to meet today's rAAV clinical requirement

(with a dosing ranging from 10^{10} to 10^{16} vg per patient, depending on the therapeutic area)

Overview of the DoE Runs



RUN1: Optimizing bioreactor operating parameters	RUN2: Optimizing cell culture parameters	RUN3: Optimizing the transfection conditions	RUN4: Optimizing the transfection conditions
DoE factors: pH <ul style="list-style-type: none"> Before TFX (Low, Middle, High) After TFX (Low, Middle, High) Stirring <ul style="list-style-type: none"> At seeding (High) Post-transfection (Low, high) Media <ul style="list-style-type: none"> Medium 1, 2 and 3 	DoE factors: Fresh Medium addition before transfection <ul style="list-style-type: none"> YES NO Media <ul style="list-style-type: none"> Medium 1 Medium 2 Medium 3 	DoE factors: Seeding VCD (10^6 cells/mL) <ul style="list-style-type: none"> Low, Middle, High DNA amount ($\mu\text{g}/10^6$ cells) <ul style="list-style-type: none"> Low, Middle, High FectoVIR[®]-AAV volume ($\mu\text{L}/10^6$ cells) <ul style="list-style-type: none"> Low, Middle, High 	DoE factors: Plasmid ratio optimization
Transfection reagent: FectoVIR [®] -AAV			
Helper plasmid: pALD-X80		Helper plasmid: pPLUS [®] -AAV Helper	
Control medium: Expi293 _{TM} Expression Medium			
		Helper plasmid control: pALD-X80	

Responses
Output / Method:
Genome Titer (vg/mL)
<ul style="list-style-type: none"> qPCR ddPCR
Total Capsid Titer (VP/mL)
<ul style="list-style-type: none"> ELISA mass-photometry
Full capsids %
<ul style="list-style-type: none"> Genome titer/Total Capsid titer mass-photometry
Transfection efficiency
efficiency (Flow cytometry)

Metabolites and ddPCR/ELISA services

Optimizing Bioreactor Operating & Cell Culture Parameters


RUN1: Optimizing bioreactor operating parameters	RUN2: Optimizing cell culture parameters
<p>DoE factors:</p> <p>pH</p> <ul style="list-style-type: none"> Before TFX (Low, Middle, High) After TFX (Low, Middle, High) <p>Stirring</p> <ul style="list-style-type: none"> At seeding (High) Post-transfection (Low, high) <p>Media</p> <ul style="list-style-type: none"> Medium 1, 2 and 3 	<p>DoE factors:</p> <p>Fresh Medium addition before transfection</p> <ul style="list-style-type: none"> YES NO <p>Media</p> <ul style="list-style-type: none"> Medium 1 Medium 2 Medium 3
Transfection reagent: FectoVIR®-AAV	
Helper plasmid: pALD-X80	Helper plasmid: pPLUS®-AAV Helper
Control medium: Expi293TM Expression Medium	
Helper plasmid control: pALD-X80	





Ambr®15

Optimizing Bioreactor Operating Parameters & Cell Culture Conditions

Run 1

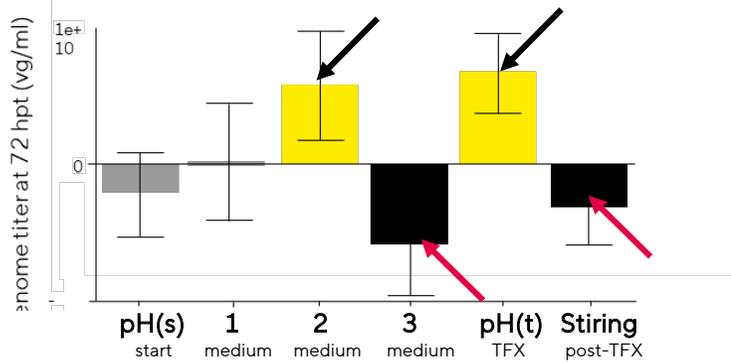


Run 2

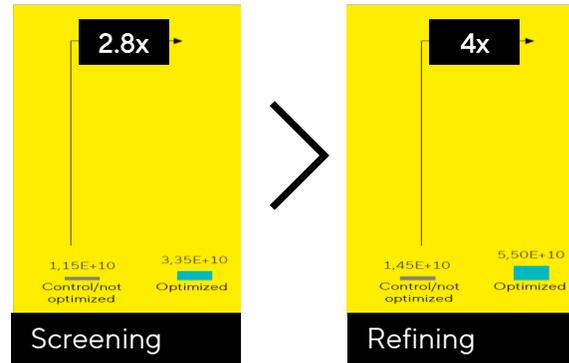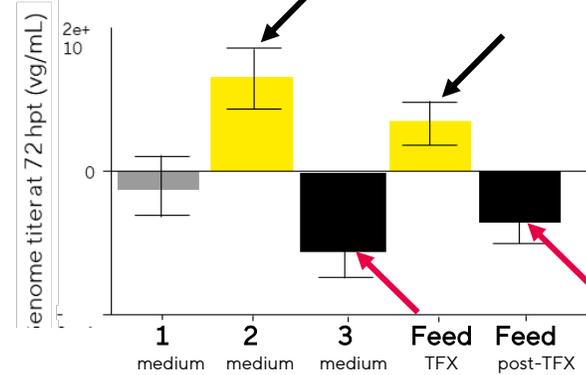

To Increase Genome Titers:

- **Medium 2** is the most performing medium.
- **Increase pH** during transfection
- **Decrease stirring speed** at the time of transfection


Optimizing Transfection conditions

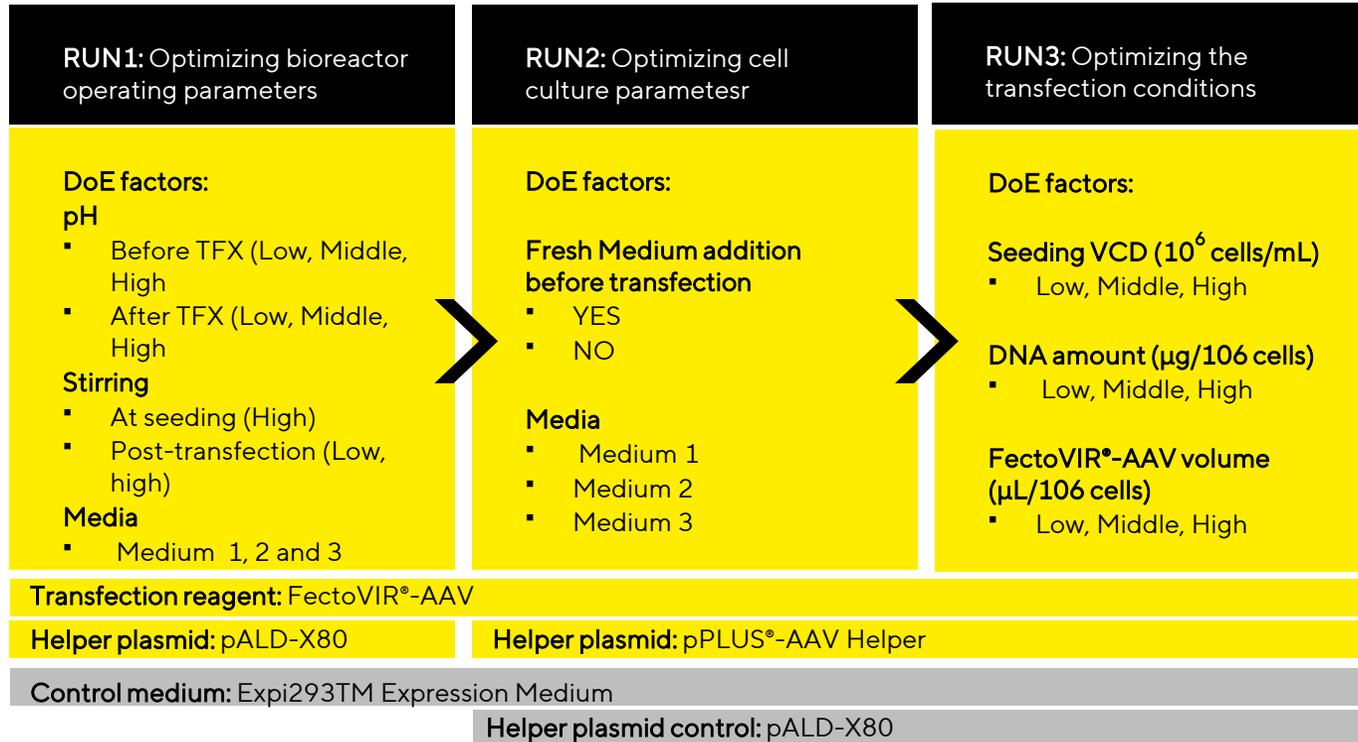

Ambr®15




Optimizing Transfection Conditions

Run 3

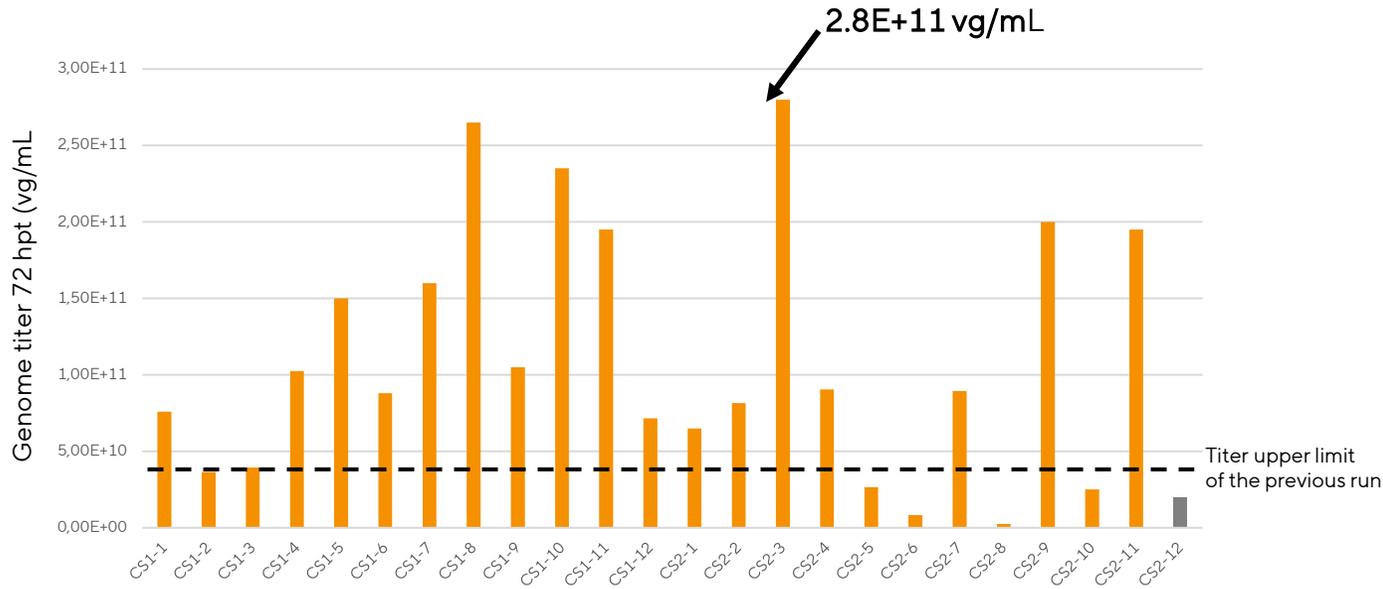


To improve productivity:

- **Decrease** cell seeding density
- **Increase** DNA amount

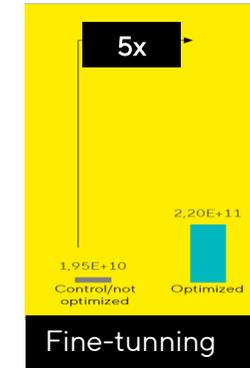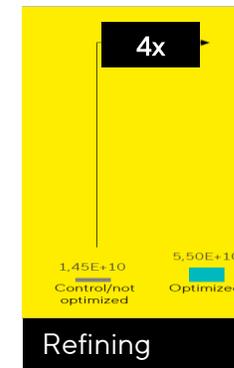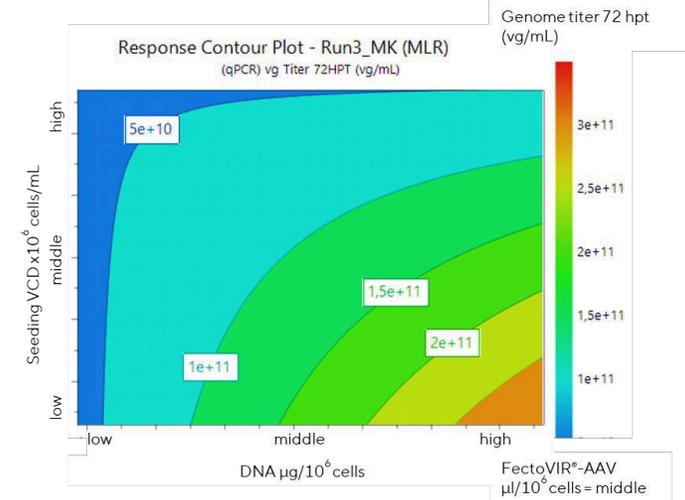


Optimizing Transfection Conditions

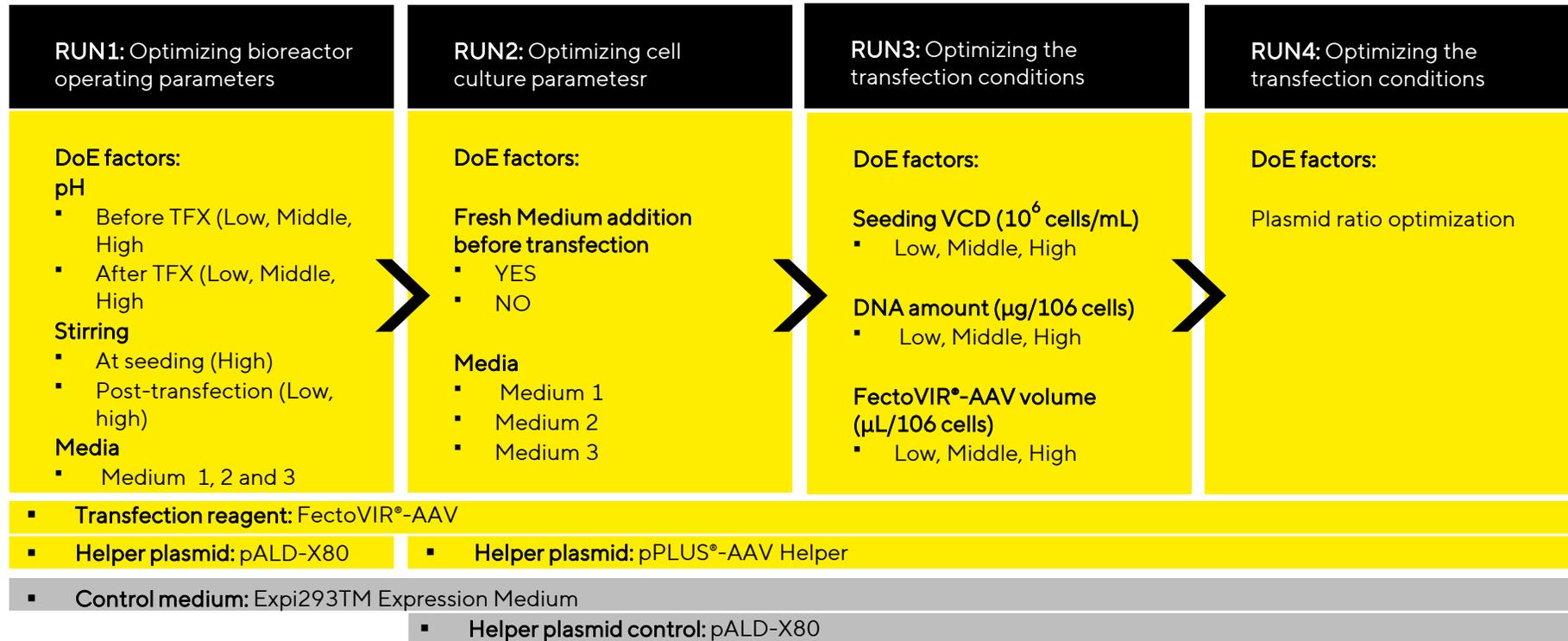




Ambr®15

Optimizing Plasmid Ratio

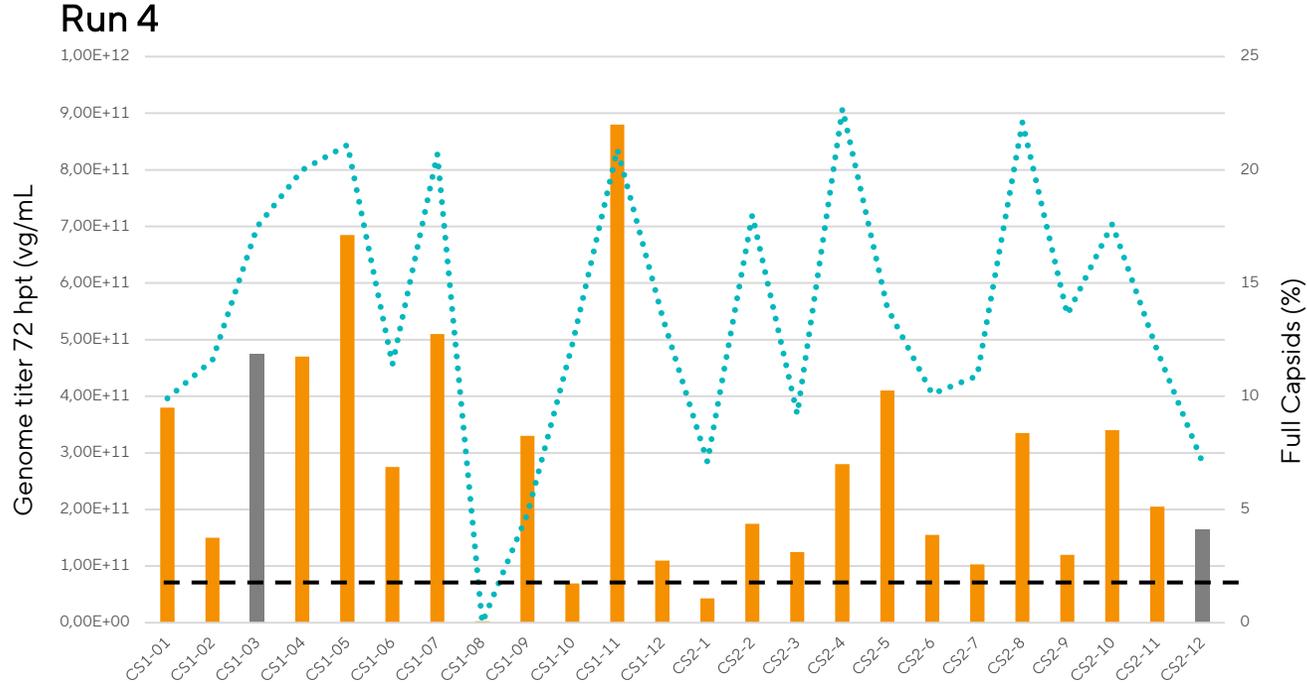


Optimal plasmid ratio for improved productivity:


- pPLUS®AAV-Helper: 45%
- pAAV-RC2: 45%
- pAAV-GFP: 10%

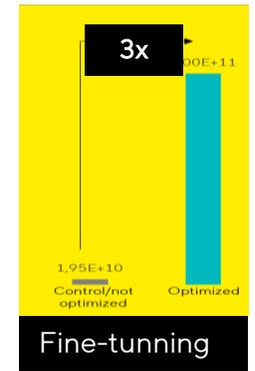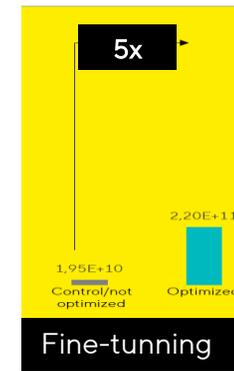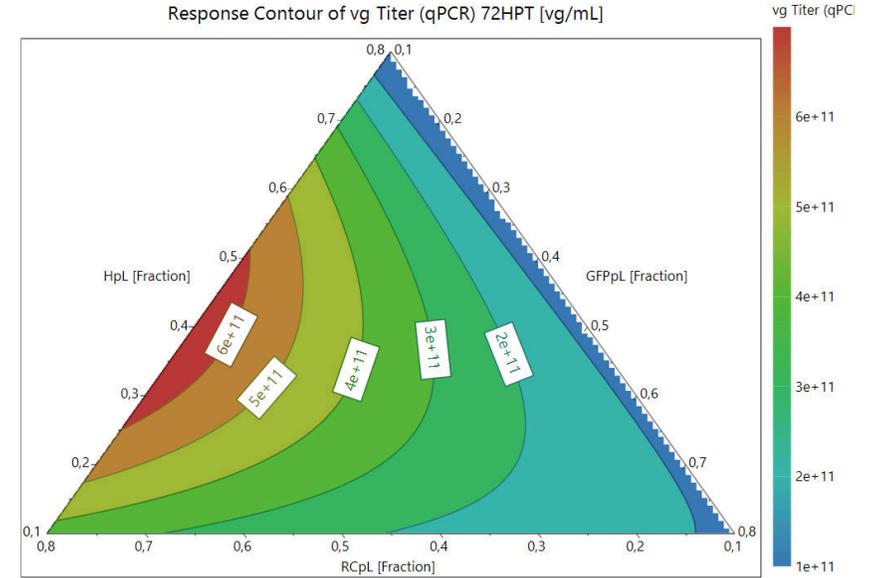

Four Experimental Designs to Achieve Significant Increase in VG

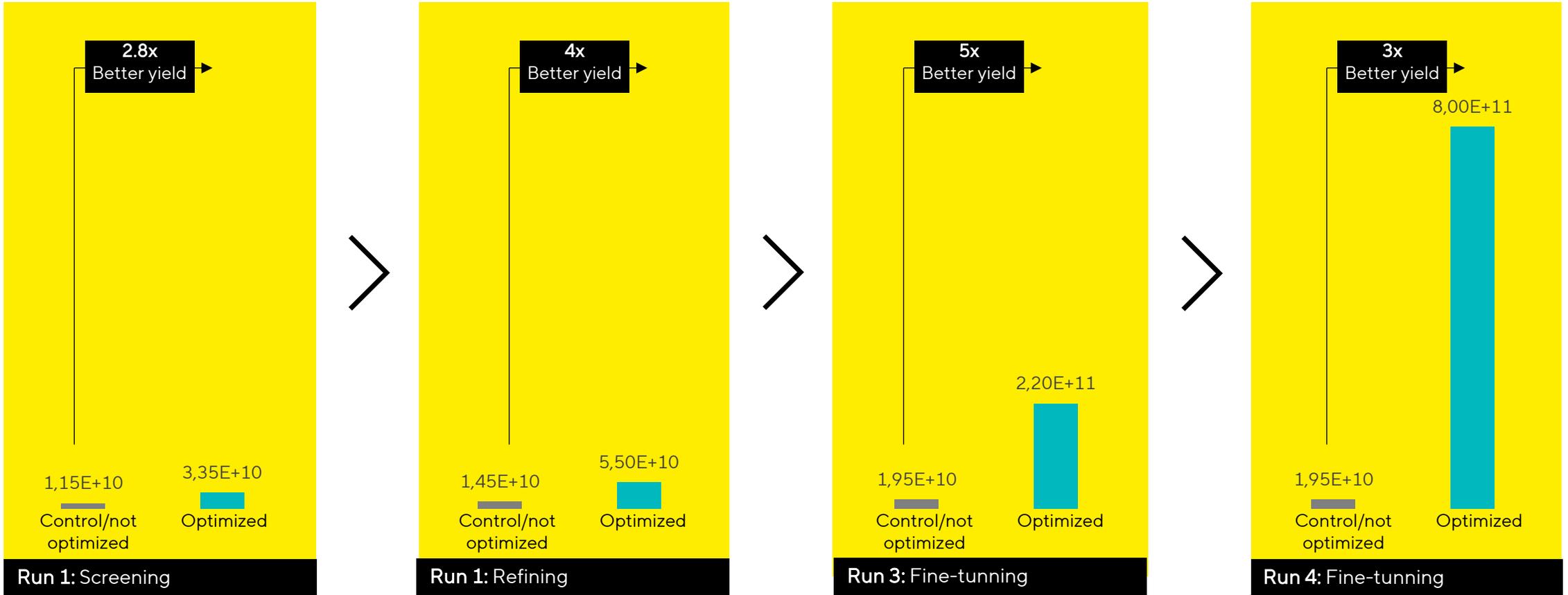



Four Experimental Designs to reach a 70-fold Increase in VG

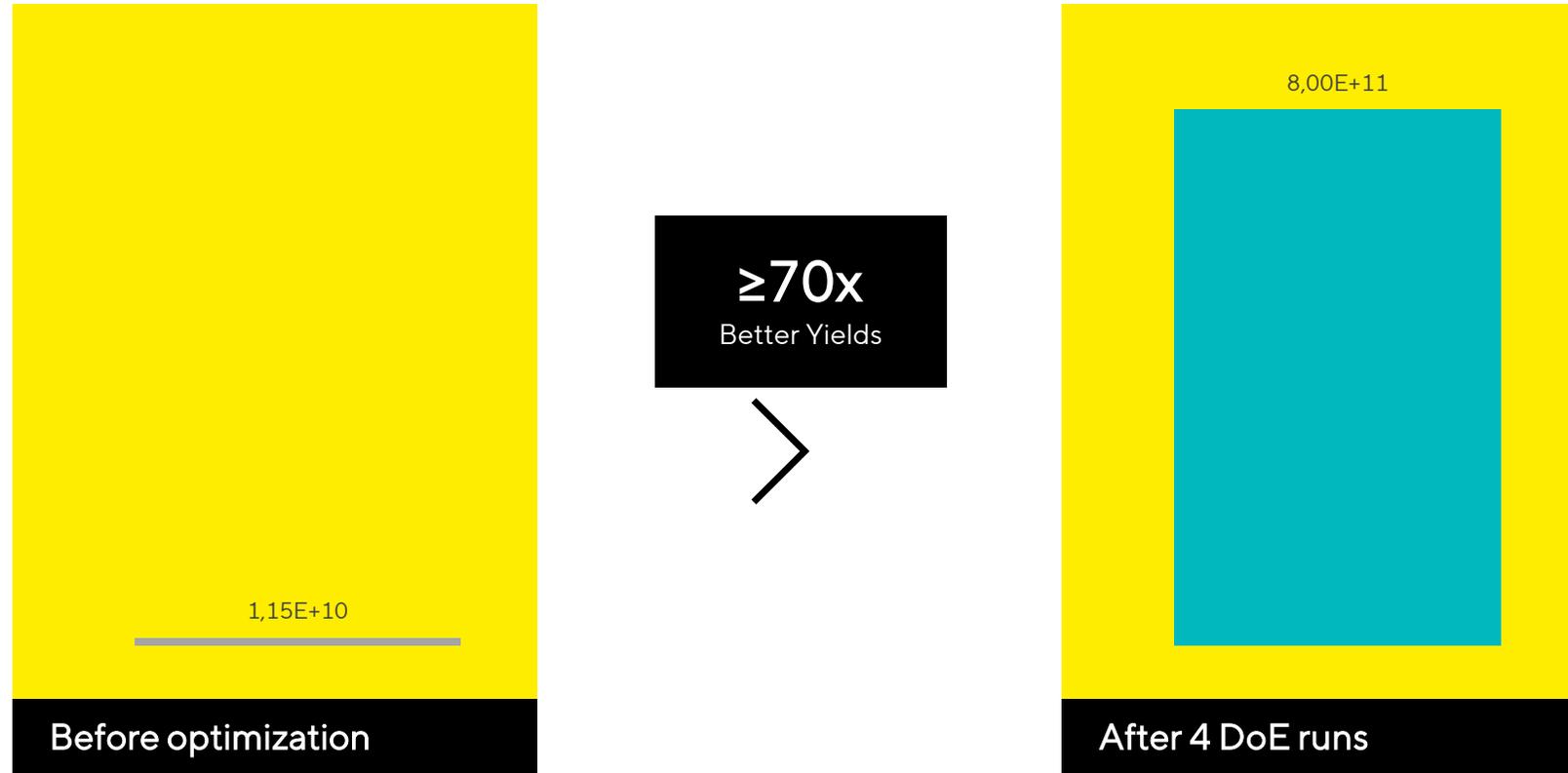

Scalable Success: Consistent Productivity at 2 L with Optimized conditions




2 L Univessel SU bioreactor

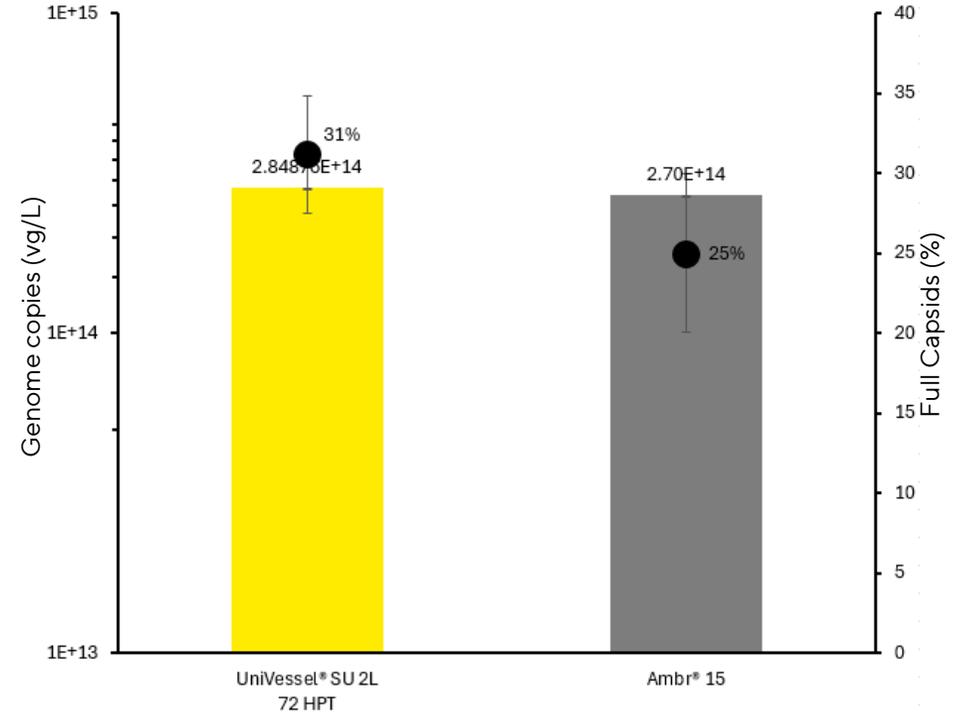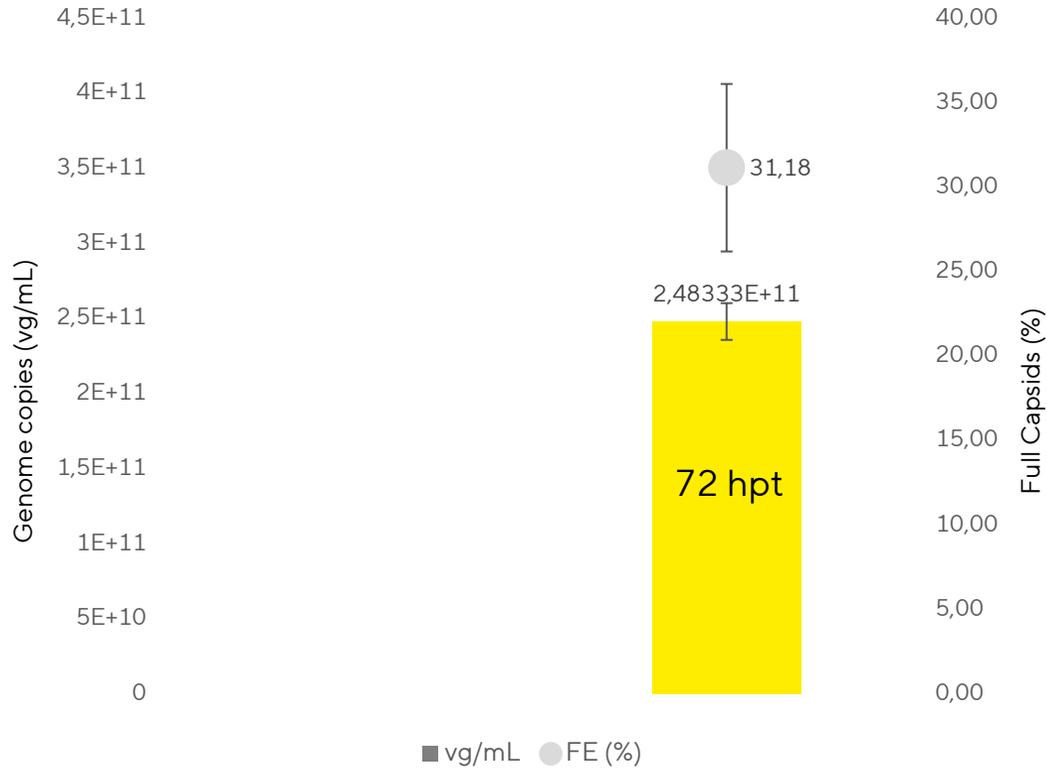

Key results and take aways

Summary and Key Takeaways

- **70-fold increase in rAAV2 productivity**
1.2E+10 → 8.8E+11 vg/mL in 4 data-driven iterations
- **Structured DoE in Ambr® 15 revealed critical multi-factor interactions**
Enabled definition of a robust operating window and deeper process understanding
- **Optimized conditions successfully translated to 2 L scale**
Demonstrating scalability and accelerated rAAV process development

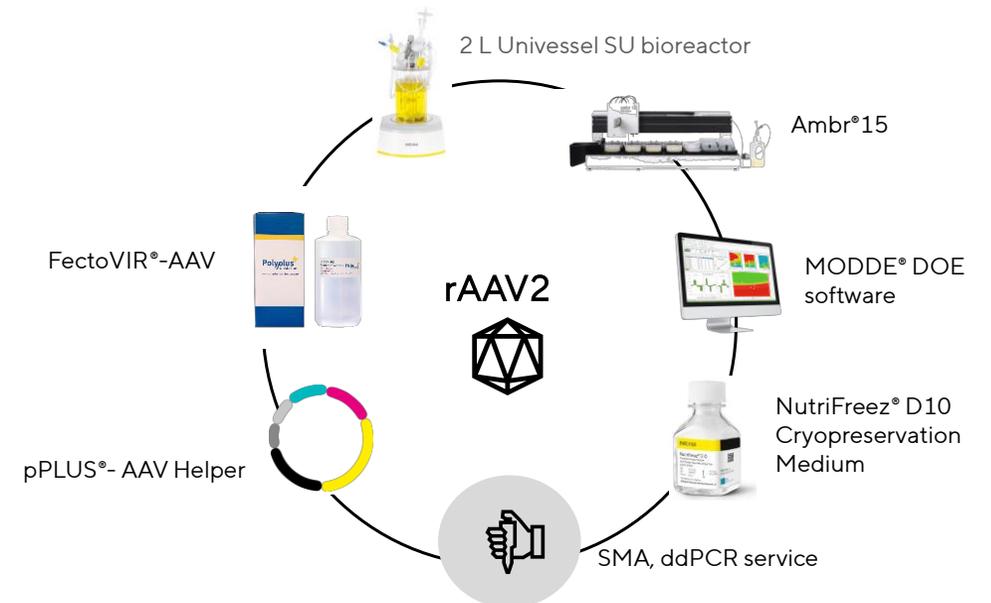

Acknowledgements



Thank You:

Érica Schulze
David Ede
Alengo Nyamay'Antu, Alengo
Maya Fuerstenau-Sharp
Geraldine Guerin-Peyrou
Noushin Dianat
Paul Giroud
Cedric Poulet
Maina Kerbrat
Thomas Jeffery
Frederique Marme-Galloo
Franziska Bollmann
Audrey LeRu
Sylvain Lebarbier
Romain Boulho

Alison Rees-Manley
Ian Ransome
Timo Schmidberger
Ryan Wray
Laure Robert
Tim Steffens
Mareike Schulz
Cedric Poulet
Mathieu Porte

Thank You:

Oumaya Adjali
Christophe Chevé
Laurence Guianvarc'h
Fabio Mirabella
Cécile Robin
Nathalie Caillé
Solène Vollant
Marie Enga
Véronique Blouin
Frédérique Broucqque
Marine Trapani
Sylvie Saleun
Emilie Audran

Booth number 79

Poster number 33

Decentralized Manufacturing of ATMPs: Navigating Regulatory Innovations and Strategic Partnerships for Enhanced Patient Care

Speaker : Ravid Grimberg

Date / Time : 18th March 2026 at 11.30-11.50 GMT

Thank you.

Specifications subject to change without notice.

© 2024 Copyright Sartorius Stedim Biotech GmbH,
August-Spindler-Strasse 11,
37079 Göttingen, Germany

Advanced Therapies 2026 | Do not replicate

The Sartorius logo is displayed in a bold, black, sans-serif font on a bright yellow rectangular background. The letters are closely spaced and have a clean, modern appearance.